

## Operational Instruction

1. **Unpack the system;**
2. **Switch the power level of the heating pad to high, and connect it to the output of the controller at back end;**
3. **Connect sensor to the controller at back end, affix sensor tip to the heating pad;**
4. **Connect the controller to electricity using the power cord provided, and turn on the controller;**
5. **Check/Set your desired temperature and duration.**

Quickly press/release SET key twice. The LED shows the current temperature setting. Use “+” and “-” keys to change the setting. When finished, press the SET again to confirm the change.

Press SET key once and release. The LED shows the current timer setting in minutes. Use “+” and “-” keys to change the setting. When finished, press the SET again to confirm the change.

6. **Wait 10-15 minutes for the heating pad temperature becomes stable.**
7. **Transfer animal onto warmed heating pad.**
8. **Insert the temperature sensor tip 1-2 cm into animal rectum for in vivo monitoring and feedback.**
9. **For more information please refer to [user's manual](#) at Docol website.**
10. **Call (888)481-0842 or email to [support@docol.com](mailto:support@docol.com) for support if you need further help.**